**[EN]: AquaVial™ Total Bacteria** detects bacteria and fungi concentrations as low as 500 CFU/ml (500 bacteria per ml) in 15 minutes at 35-40 °C (95-104 °F), or in 30 minutes at 20-25 °C (68-77 °F).

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**[EN]: IMPORTANT:**

Please read the directions carefully and watch the instructional video, before starting the test.

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**[EN]: DISPOSAL:**

T-Cup components, are recyclable. Please recycle whenever possible. Vials can be disposed of in the trash.

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**[EN] DIRECTIONS FOR USE:**

**VIDEO INSTRUCTIONS:** [www.aquavial.ca](http://www.aquavial.ca)


**Step 1: Collect and Concentrate Water Sample**

**OPEN T-CUP:** Separate the top and bottom halves of the T-Cup by twisting the top half counter clockwise.

**ADD WATER SAMPLE:** Fill the bottom of the T-Cup halfway with the sample water.

**CONCENTRATE:** Twist top of the T-Cup clockwise into bottom section until it stops, forcing water through the integrated T-Cup filter. Water will pour through and spill over the sides of the plunger. This is normal. The T-Cup is intended to sample this exact amount of water.

**DISCARD:** Drain excess water from the top and sides of the T-Cup.

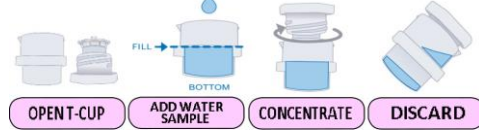**Step 2: Add Distilled Water**

This step is required to ensure the water sample collected and concentrated in Step 1 is washed of chemicals, like chlorine, used to treat the water, which may affect your test results. The addition of distilled water will not dilute or interfere with the water sample to be tested. The T-Cup system has a built-in filter that retains the bacteria at the bottom, while eliminating chemicals through the filter.

**OPEN T-CUP:** Open the T-Cup by twisting the top half counter-clockwise.

**ADD DISTILLED WATER:** Carefully tear the top section of the distilled water pouch and pour distilled water into the bottom half of the T-Cup.

**CONCENTRATE:** Twist the top of the T-Cup into the bottom clockwise to close the T-Cup.

**DISCARD:** Pour out any excess water from the top of the T-Cup.

**Step 3: Test Water Sample**


**OPEN T-CUP:** Open the T-Cup by twisting the top half counter-clockwise.

**SNAP OPEN VIAL:** Hold AquaVial™ Total Bacteria vial (pink solution) into your hand firmly in a vertical position with the tip of the cap just above your index finger. Apply side pressure with your thumb to snap the cap open. **BE CAREFUL NOT TO TILT OR SPILL AQUAVIAL™**

**FILL AQUAVIAL™:** Pour the unfiltered water from the bottom of the T-Cup into the AquaVial™ Total Bacteria vial. In most cases the liquid will not reach the fill line marked on the vial, which is normal. Sample water should not exceed the fill line; or stop filling water in the vial after this line. Close AquaVial™ Total Bacteria vial using one of the two plug caps provided.


**Tip:** Take a picture of the vial against a white background.

**INCUBATE:** Let the AquaVial™ Total Bacteria vial incubate in an upright position for 30 minutes and check for changes in colour. If you are not sure if the vial changed color, let the vial incubate another 30 minutes and check color again.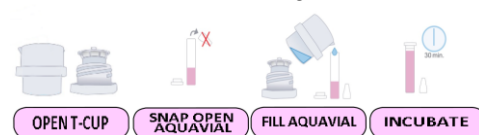

Compare the color of the reagent in the test vial against the color diagram below:

35-40°C (95-104°F) – 15 min OR 20-25°C (68-77°F) – 30 min

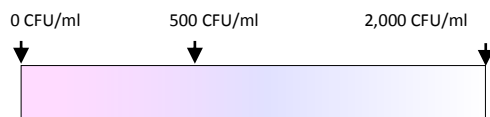If the solution changes its colour to clear, **extreme caution should be exercised**, and a sample should be taken to a laboratory for further testing. This test is intended for **rapid screening only**.

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